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**Mascot Database Manager Configuration Import**

**Note:**  
Please disable other programs and scripts that configure or update sequence databases before continuing.

Database Manager can keep your sequence database configuration up to date and schedule the download of new files.

If you choose to use Database Manager, you should not edit mascot.dat in future, because your changes will be lost. If you do not wish to use Database Manager now, please [return to the Configuration Editor](#).

Database Manager has two kinds of database definitions:

1. *Predefined database definition*: Automatic configuration by downloading up-to-date information from the Matrix Science website.
2. *Custom database definition*: Manually maintained configuration.

You can choose to keep the existing definition (Keep as Custom) or synchronise with a predefined definition (Synchronise). Synchronisation means the predefined database definition will be applied, and the database files (FASTA, reference, taxonomy) will be updated.

If you do not want to update the files now, choose Keep as Custom. You can synchronise at any time in the future by choosing Synchronise Custom Definitions from the Database menu.

Database definition	Action	Match quality
SwissProt	<input type="radio"/> Keep as Custom <input checked="" type="radio"/> Synchronise with SwissProt_ID <input type="radio"/> Synchronise with SwissProt_AC poor	perfect Differences

Import database configuration by choosing Import.

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proxy settings」を

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**Proxy settings**

You may need to define proxy settings to enable Database Manager to access the Internet, depending on your local network and Internet connection.

HTTP proxy strategy:

1. Try Database Manager proxy settings (currently no proxy settings defined).
2. Try mascot.dat settings (currently no proxy settings defined).
3. Try proxy settings defined in the system environment (currently no proxy settings defined for CGI scripts).
4. Direct connection without proxy.

URL test **succeeds**; Database Manager can access [http://www.matrixscience.com/downloads/databases\\_1.xml](http://www.matrixscience.com/downloads/databases_1.xml) from the Mascot server ma

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### Sequence directory settings

Sequence databases are usually stored in a single sequence directory. The default path is `C:/inetpub/mascot/sequence`, but this can be changed below.

Note that you can always change the sequence directory for individual database definitions when creating a new database definition.

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(If you leave this field blank, the sequence directory will be reset to the default path.)

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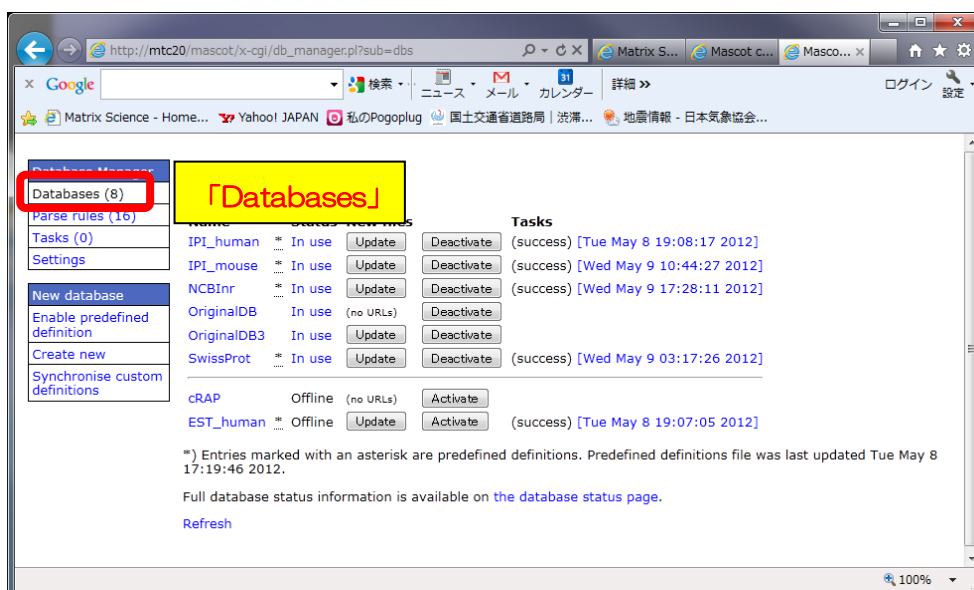
**Databases**

Name	Status	New files	Tasks
IPI_human	* In use	Update	Deactivate (success) [Tue May 8 19:08:17 2012]
IPI_mouse	* In use	Update	Deactivate (success) [Wed May 9 10:44:27 2012]
NCBINr	* In use	Update	Deactivate (success) [Wed May 9 17:28:11 2012]
OriginalDB	In use	(no URLs)	Deactivate
OriginalDB3	In use	Update	Deactivate
SwissProt	* In use	Update	Deactivate (success) [Wed May 9 10:44:27 2012]
cRAP	Offline	(no URLs)	Activate
EST_human	* Offline	Update	Activate (success) [Tue May 8 19:07:05 2012]

\*) Entries marked with an asterisk are prepared by MASCOT. Full database status information is available in the database file.

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Matrix Science - Home x Mascot configuration x Mascot Database Ma... x

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**Most recent finished task**  
[Wed May 9 03:17:26 2012] (success) No new FASTA file available; database SwissProt is up to date.

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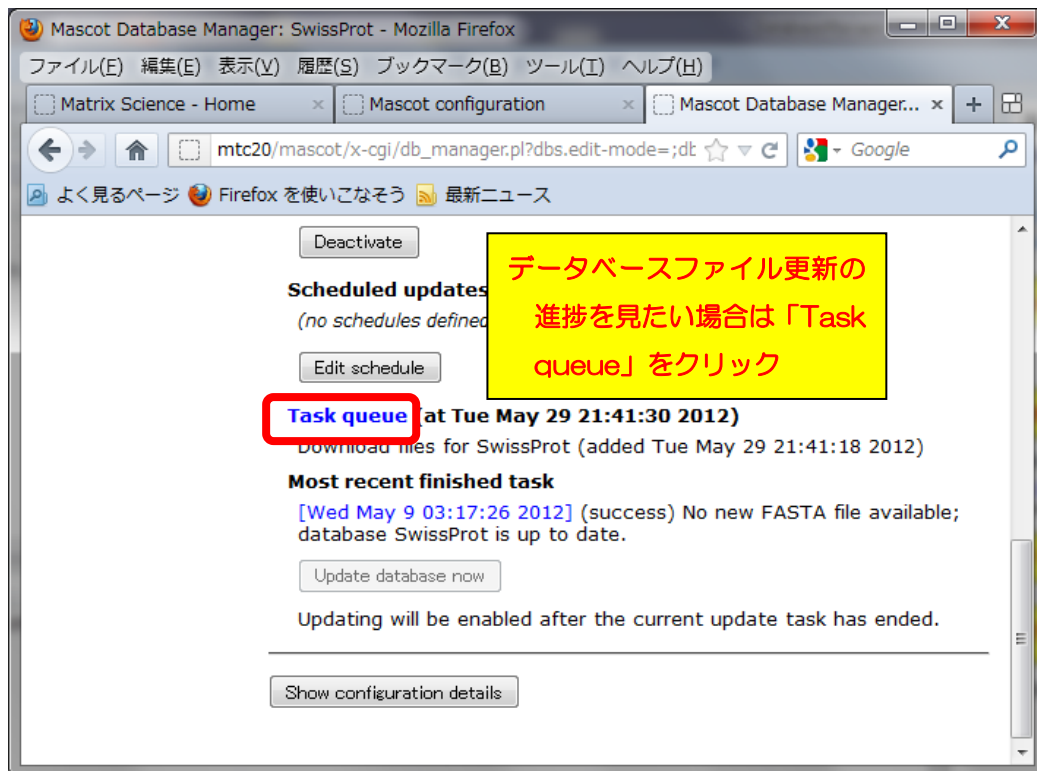
① 「databases」

Databases

Name	Status	New files	Tasks
IPI_human	In use	Update Deactivate	(success) [Tue May 8 19:08:17 2012]
IPI_mouse	In use	Update Deactivate	(success) [Wed May 9 10:44:27 2012]
NCBINr	In use	Update Deactivate	(success) [Wed May 9 17:28:11 2012]
OriginalDB	In use	Update Deactivate	(success) [Wed May 9 03:17:26 2012]
SwissProt	In use	Update Deactivate	(success) [Wed May 9 03:17:26 2012]

② database 選択

Deactivate

Scheduled updates  
(no schedules defined)

Edit schedule

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Task queue (at Tue May 29 21:41:30 2012)  
Download files for SwissProt (added Tue May 29 21:41:18 2012)

Most recent finished task  
[Wed May 9 03:17:26 2012] (success) No new FASTA file available; database SwissProt is up to date.

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**Filename pattern**  
Uniprot\_Human\_\*.fasta

**No local FASTA file**

Database files must be present before the database can be configured and activated. You have the following options:

- Download from remote URL or...
- Upload file using web browser
- Copy file manually

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Edit filename pattern

Please,

- copy the FASTA file to C:/inetpub/mascot/sequence/Uniprot\_Human/current;
- rename the file to match the filename pattern, e.g. Uniprot\_Human\_20150512.fasta (in the pattern, \* matches any string of characters);
- and refresh this page to continue.

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The image shows two screenshots of the Mascot database manager interface. The top screenshot displays the configuration page for the 'Uniprot\_Human' database. The 'Database status' section shows the database is 'Offline' and an 'Activate' button is highlighted with a red circle. A yellow callout box points to this button with the text: ファイル名が認識されたら「Activate」.

The bottom screenshot shows the 'Database Status' page. It lists the status of several databases. The 'Uniprot\_Human' database is highlighted with a red box, and a yellow callout box points to it with the text: Database Status 画面でデータベース構築確認.

**Database: Uniprot\_Human**

**Database Manager**

- Databases (12)
- Parse rules (21)
- Tasks (0)
- Settings

**Uniprot\_Human**

- Copy
- Delete

**Name**  
Uniprot\_Human

**Database type**  
Amino acid (protein database)

**Database directory**  
C:/inetpub/mascot/sequence/Uniprot\_Human/current

**Filename pattern**  
Uniprot\_Human\_\*.fasta

**Matching files in directory**  
Uniprot\_Human\_20120530.fasta (43.72 MB)

**Database status**  
Offline

**Activate**

**Scheduled updates**  
(no schedules defined)

**Database Status**

Filename = Environmental\_ES1\_111.fasta Pathname = C:/inetpub/mascot/sequence/Environmental\_ES1/Current  
Status = In use [Statistics](#) [Unidentified taxonomy](#)  
State Time = Wed May 30 11:31:57 # searches = 0  
Mem mapped = YES Request to mem map = YES Request unmap = NO Mem locked = NO  
Number of threads = -1 Current = YES

Name = IPI\_bovine Pathname = C:/inetpub/mascot/sequence/IPI\_bovine/Current  
Filename = IPI\_bovine\_3.73.fasta Pathname = C:/inetpub/mascot/sequence/IPI\_bovine/Current  
Status = In use [Statistics](#)  
State Time = Wed May 30 11:32:16 # searches = 0  
Mem mapped = YES Request to mem map = YES Request unmap = NO Mem locked = NO  
Number of threads = -1 Current = YES

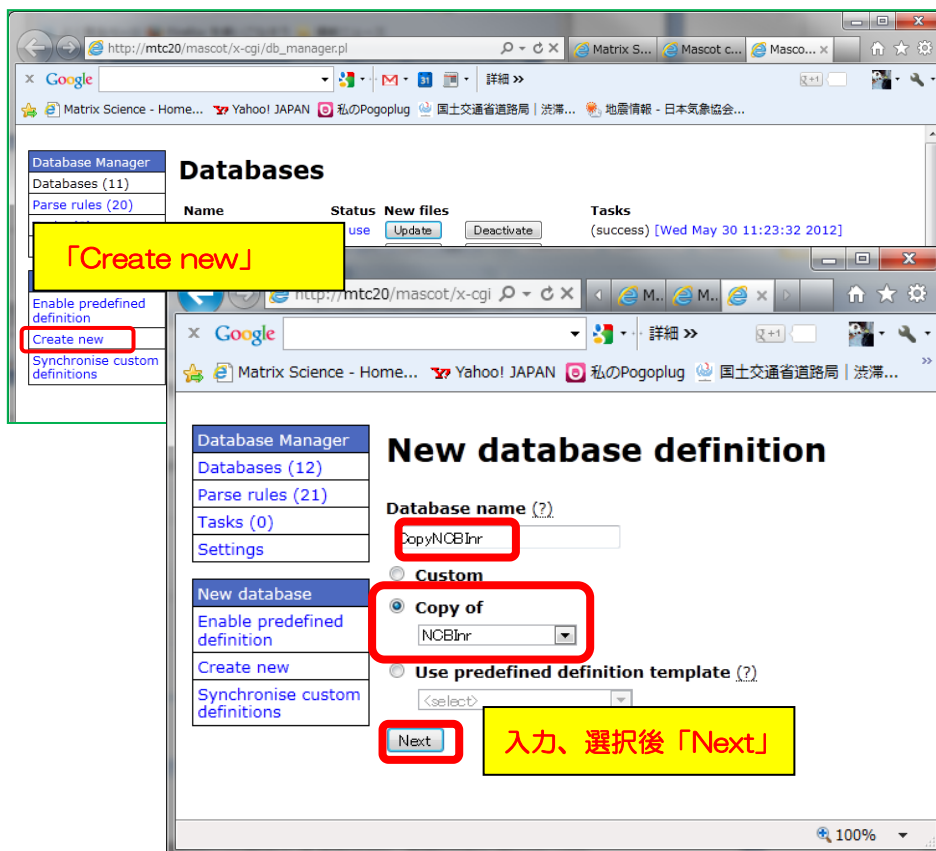
Name = Uniprot\_Human Pathname = C:/inetpub/mascot/sequence/Uniprot\_Human/current/Uniprot\_Human\_20120530.fasta  
Filename = Uniprot\_Human\_20120530.fasta Pathname = C:/inetpub/mascot/sequence/Uniprot\_Human/current/Uniprot\_Human\_20120530.fasta  
Status = Creating compressed files 39% complete  
State Time = Wed May 30 14:15:40 # searches = 0  
Mem mapped = NO Request to mem map = YES Request unmap = NO Mem locked = NO  
Number of threads = -1 Current = NO

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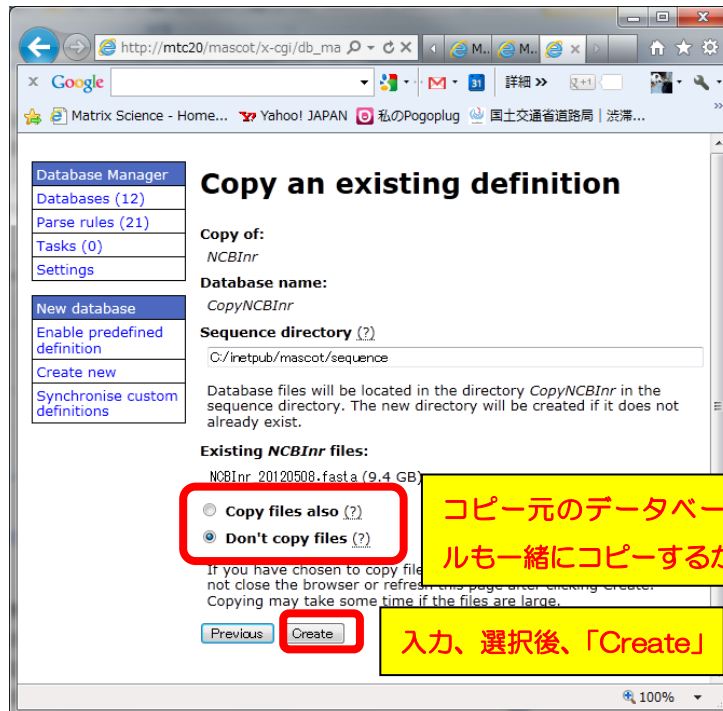
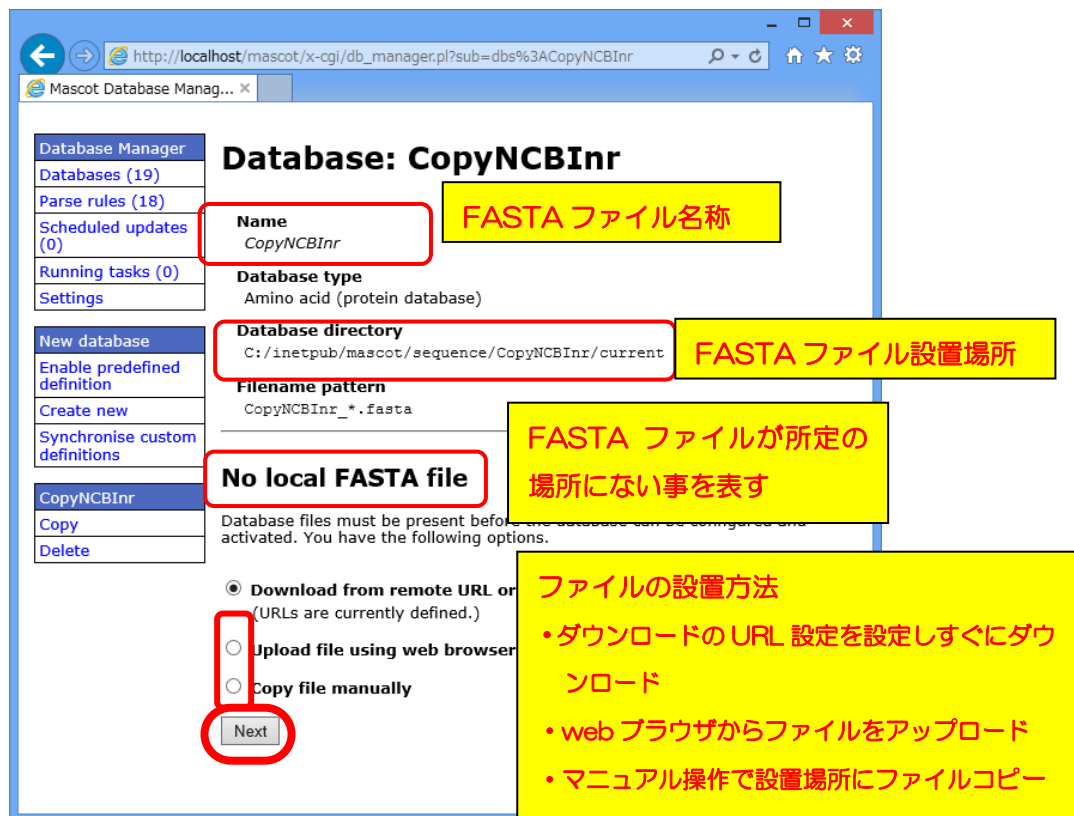
The screenshot shows the Mascot Database Manager interface. The main content area displays the following information for the 'CopyNCBIInr' database:

- Name:** CopyNCBIInr
- Database type:** Amino acid (protein database)
- Database directory:** C:/inetpub/mascot/sequence/CopyNCBIInr/current
- Filename pattern:** CopyNCBIInr\_\*.fasta

The 'Copy files manually' section includes the following instructions:

- copy the FASTA file to C:/inetpub/mascot/sequence/CopyNCBIInr/current;
- rename the file to match the filename pattern, e.g. CopyNCBIInr\_20150512.fasta in the pattern. \* matches any string of characters);
- and refresh this page to contin

Three yellow callout boxes with red text provide additional instructions:

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
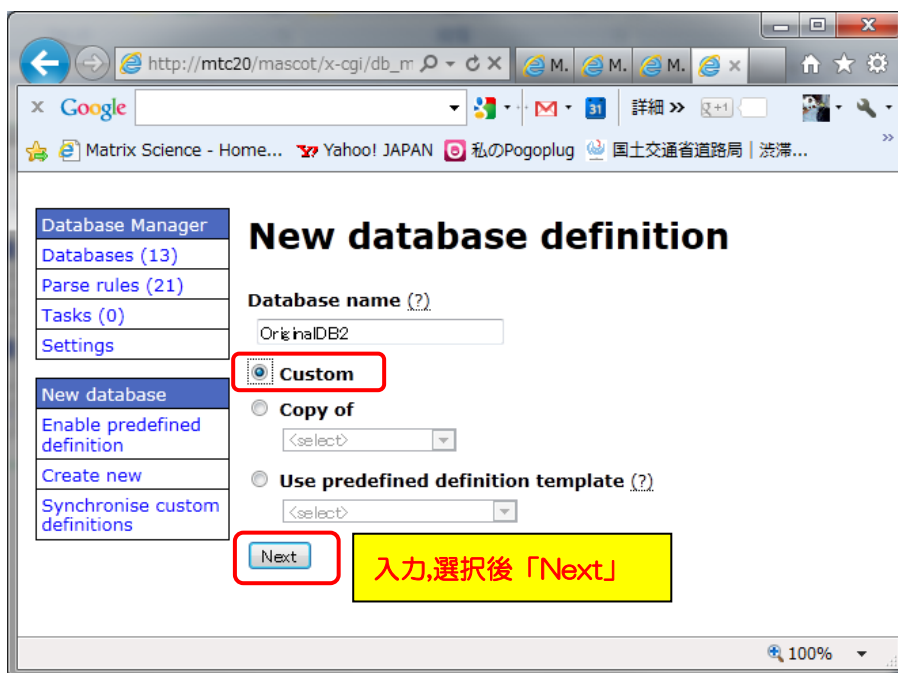
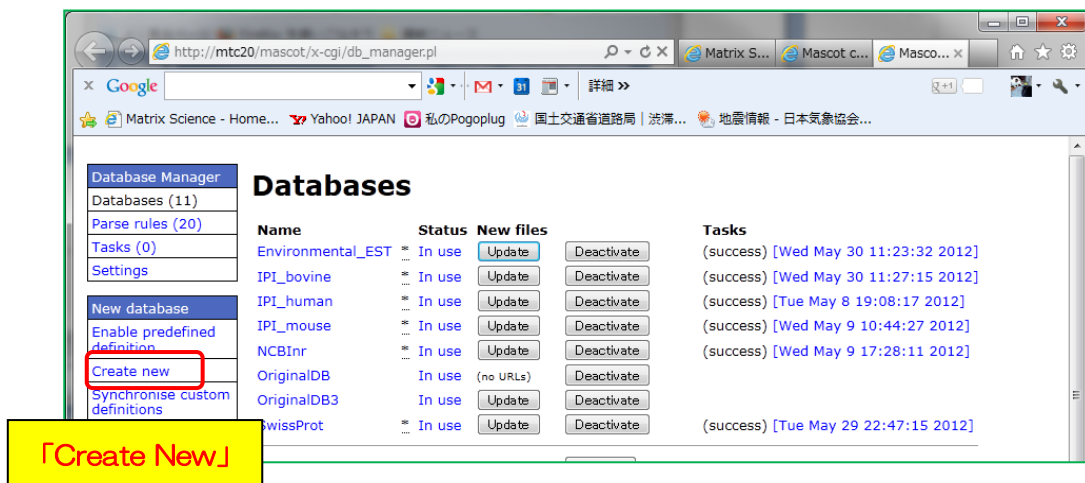
The screenshot shows the 'Custom definition' page in the Mascot Database Manager. The page is titled 'Custom definition' and has a sidebar on the left with navigation links: Database Manager, Databases (16), Parse rules (17), Scheduled updates (0), Running tasks (0), Settings, New database, Enable predefined definition, Create new, and Synchronise custom definitions. The main content area includes the following fields and options:

- Database name:** OriginalDB2
- Database type:** Radio buttons for 'Amino acid (protein database)' (selected) and 'Nucleic acid (DNA database)'. An annotation points to this section: 「Database type」 アミノ酸か核酸か
- Sequence directory (?):** Text input field containing 'C:/inetpub/mascot/sequence'. An annotation points to this field: 「Sequence directory」 配列データベースファイルの置き場所
- Download or copy files automatically (?):** Radio button (selected) for automatic download. Below it are three text input fields for 'FASTA file URL or path to source file on Mascot Server hard disk', 'Version file URL or path to source file on Mascot Server hard disk', and 'Reference file URL or path to source file on Mascot Server hard disk'. Each has a checkbox for 'Delete original file after copying'. An annotation points to the 'Download database files' option: 「Download database files」 自動更新プログラムにおけるファイルの取得先
- Upload or copy files manually (?):** Radio button (unselected). An annotation points to this option: 「Upload or Copy files manually」 上記設定にて自動ダウンロードしない場合に選択

At the bottom of the page, there are two buttons: 'Previous' and 'Create'. The 'Create' button is highlighted with a red box and an annotation: 入力、選択後「Create」

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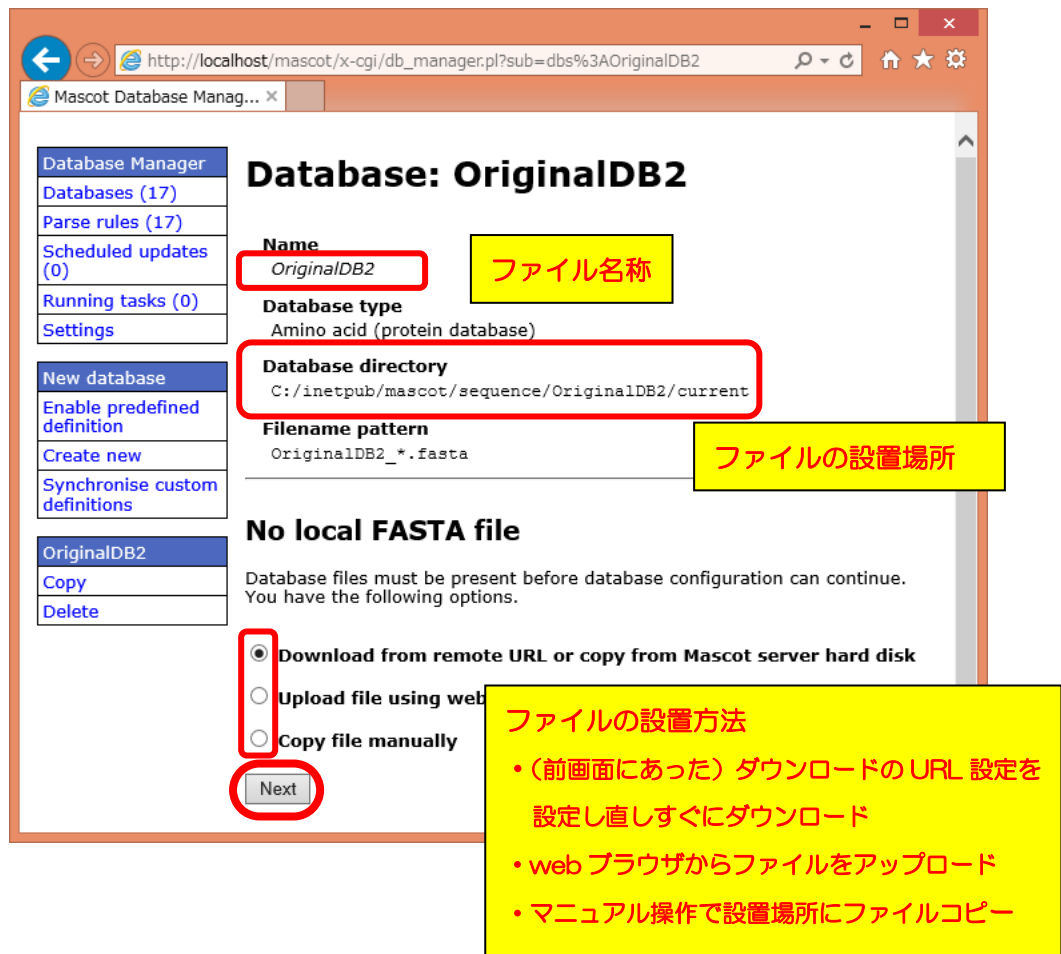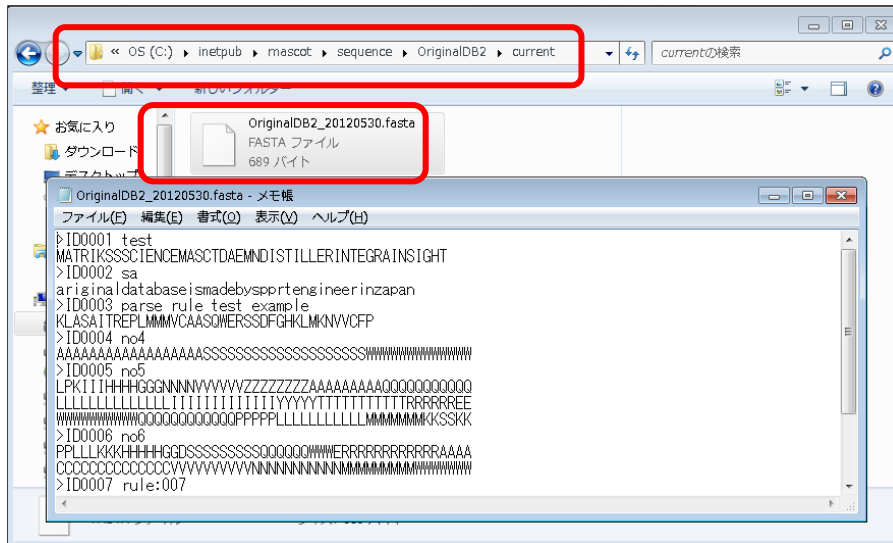
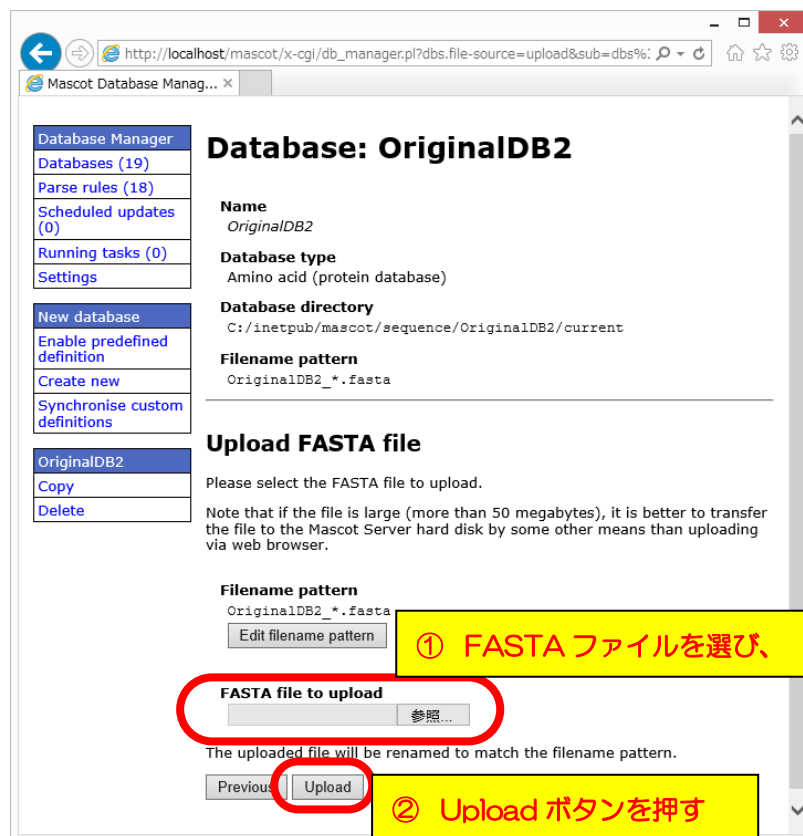
**Choose** ① Accession の「Choose」

**Description parse rule (?)**  
(none)

**FASTA file accession parse rule for OriginalDB2**

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**Parse rules (5 suitable)**

Parse rule	Match	Extracted data
<input type="radio"/> >[^ ]* ¥(.*)	5/5	1. test ▶ 4 more matches
<input type="radio"/> >¥(.*)	5/5	1. ID0001 test ▶ 4 more matches
<input checked="" type="radio"/> >¥([^\ ]*)	5/5	1. ID0001 ▶ 4 more matches
<input type="radio"/> >¥([^\ ]*)	5/5	1. ID0001 test ▶ 4 more matches
<input type="radio"/> >¥([^\ ]*)	5/5	1. >ID0001 test ▶ 4 more matches

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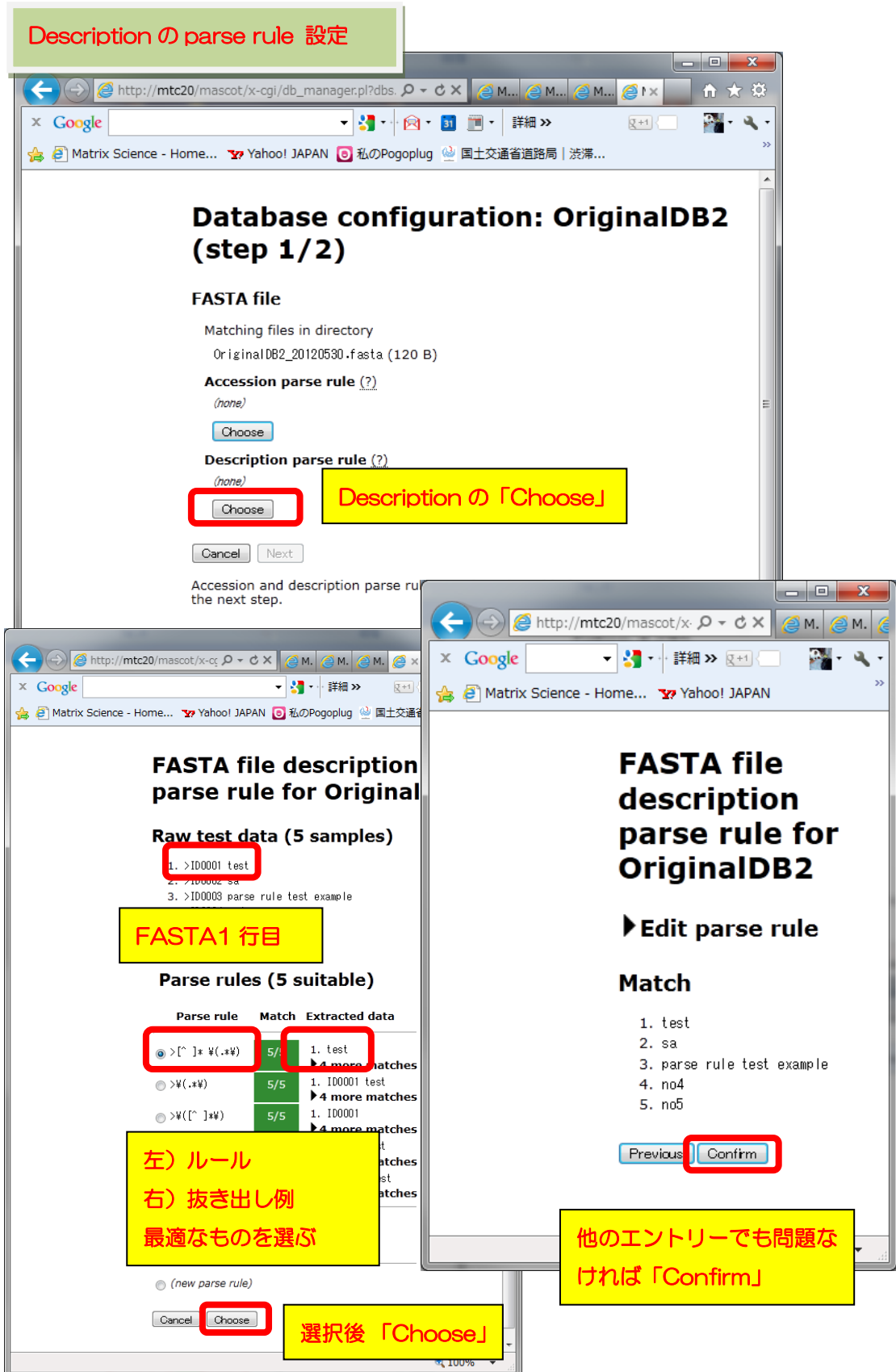
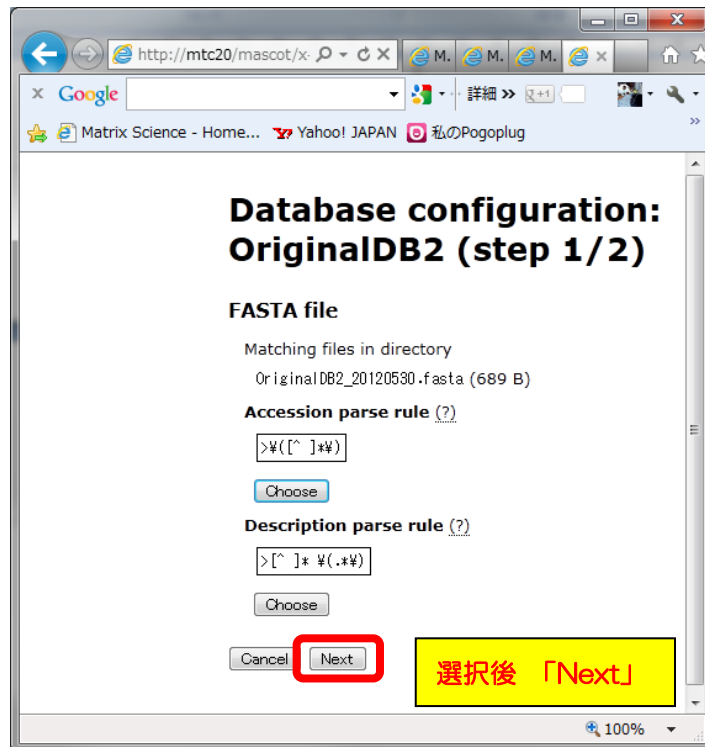
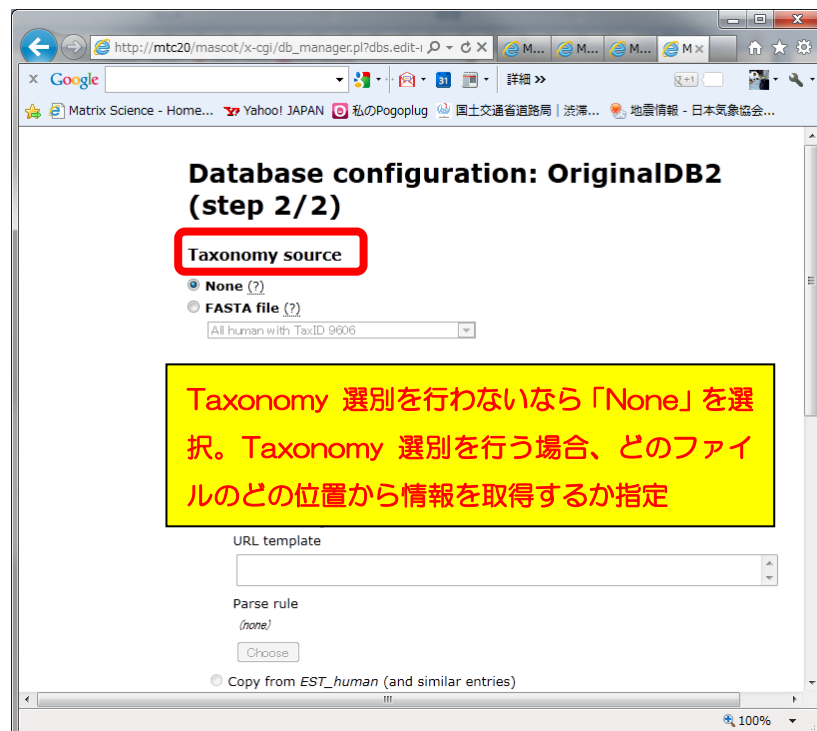
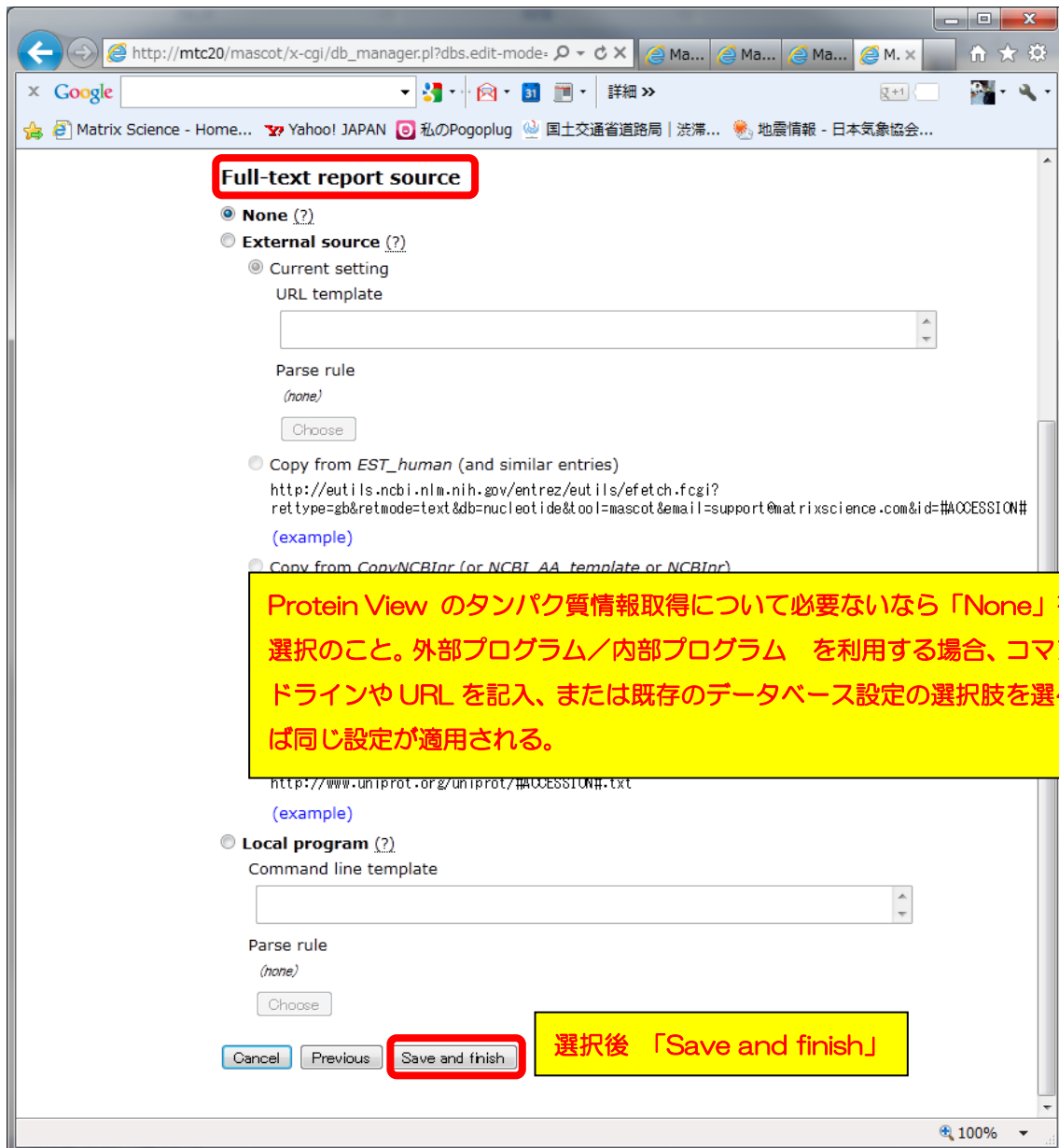
The image shows two screenshots of the Mascot database manager interface. The top screenshot displays the configuration page for 'Database: OriginalDB2'. The 'Database status' section shows 'Offline' and an 'Activate' button, which is highlighted with a red box. A yellow callout box next to it says '「Activate」ボタンで使用開始'. Below this, the 'Scheduled updates' section shows '(no schedules defined)' and an 'Update database now' button. The bottom screenshot shows the 'ms-stat' status page. It lists the status of various databases. The 'OriginalDB2' entry is highlighted with a red box and has a yellow callout box that says 'Database 構築開始'. The status for 'OriginalDB2' is 'Creating compressed files 97% complete'.

**Database: OriginalDB2**

**Name**  
OriginalDB2

**Database type**  
Amino acid (protein database)

**Database directory**  
C:/inetpub/mascot/sequence/OriginalDB2/current

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FASTA file URL has not been defined; database updates are disabled.

Mem mapped = YES Request to mem map = YES Request unmap = NO Mem locked = NO  
Number of threads = -1 Current = YES

Name = Uniprot\_Human Family = C:/inetpub/mascot/sequence/Uniprot\_Human/current  
Filename = Uniprot\_Human\_20120530.fasta Pathname = C:/inetpub/mascot/sequence/Uniprot\_Humar  
Status = In use [Statistics](#)

State Time = Wed May 30 14:15:52 # searches = 0  
Mem mapped = YES Request to mem map = YES Request unmap = NO Mem locked = NO  
Number of threads = -1 Current = YES

**Database 構築開始**

Name = OriginalDB2 Family = C:/inetpub/mascot/sequence/OriginalDB2/current/  
Filename = OriginalDB2\_20120530.fasta Pathname = C:/inetpub/mascot/sequence/OriginalDB2/cur  
Status = Creating compressed files 97% complete  
State Time = Wed May 30 17:13:58 # searches = 0  
Mem mapped = NO Request to mem map = YES Request unmap = NO Mem locked = NO  
Number of threads = -1 Current = NO

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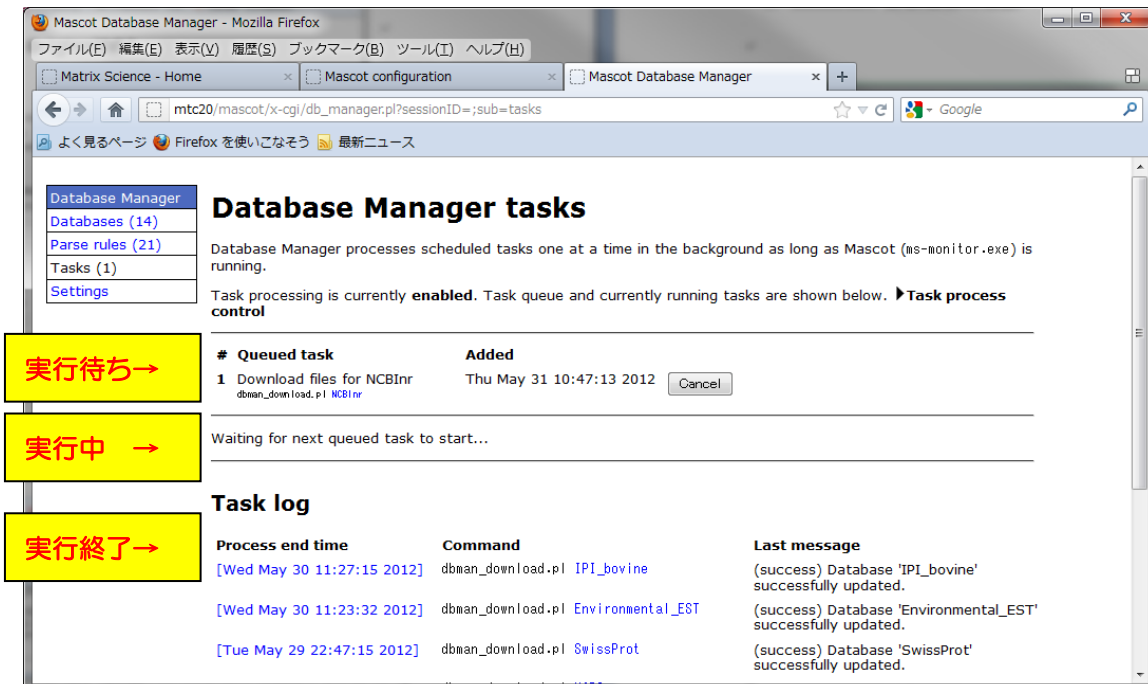
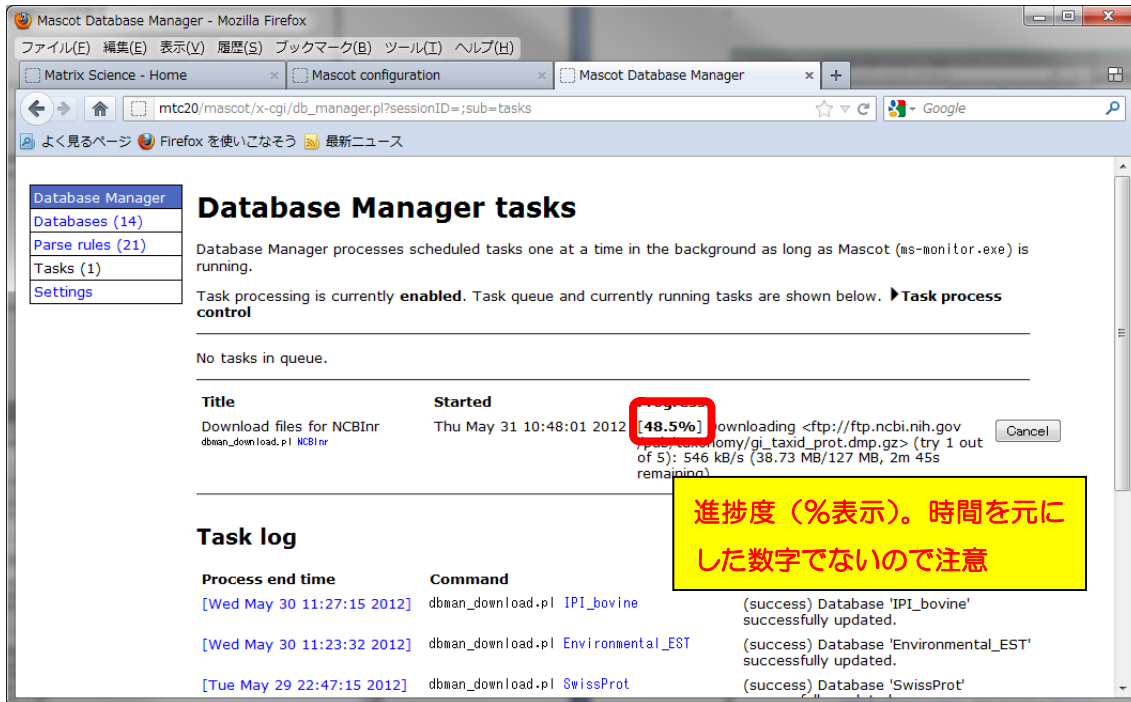
The screenshot shows the Mascot Database Manager interface in a Mozilla Firefox browser. The page title is "Database Manager tasks". The main content area displays a task in progress: "Download files for NCBItr" with a progress bar at 48.5%. The task details include the start time "Thu May 31 10:48:01 2012" and the command "dbman\_download.pl NCBItr". A yellow callout box highlights the progress percentage with the text: "進捗度 (%表示)。時間を元にした数字でないので注意". Below the task list is a "Task log" section showing previous successful updates for databases "IPI\_bovine", "Environmental\_EST", and "SwissProt".

Title	Started	Progress	Details
Download files for NCBItr <small>dbman_download.pl NCBItr</small>	Thu May 31 10:48:01 2012	48.5%	downloading <ftp://ftp.ncbi.nih.gov/pub/blast/blast_taxonomy/gi_taxid_prot.dmp.gz> (try 1 out of 5): 546 kB/s (38.73 MB/127 MB, 2m 45s remaining)

Process end time	Command	Details
[Wed May 30 11:27:15 2012]	dbman_download.pl IPI_bovine	(success) Database 'IPI_bovine' successfully updated.
[Wed May 30 11:23:32 2012]	dbman_download.pl Environmental_EST	(success) Database 'Environmental_EST' successfully updated.
[Tue May 29 22:47:15 2012]	dbman_download.pl SwissProt	(success) Database 'SwissProt' successfully updated.

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